



A colossal CONSERVATION EXPERIENCE

Name: _____

INGREDIENTS



Pup's Saliva



Dish Soap



Salt



Rubbing Alcohol



Toothpick



Clear Dish

INSTRUCTIONS FOR EXTRACTING THE PUP DROOL

1



Gently collect the drool from the dog's lips
(but you need to have a drooly dog!)

2



Put the drool in a clear dish.

3




Add an equal amount of dish soap
to the drool.

4



Add a pinch of salt.

5

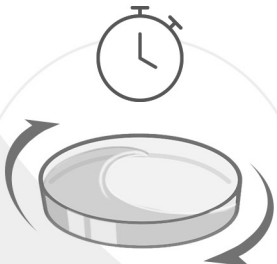

Gently swirl to mix and let sit for
a minute.

6



Gently pour rubbing alcohol along
the side of dish.

7



Watch the white, stringy strands
of DNA forming.

8



Use a toothpick to lift the sticky
strands of DNA.

